


# Frequency Dependence of Dielectric Constant

SES Instruments Pvt Ltd.

FDD-01

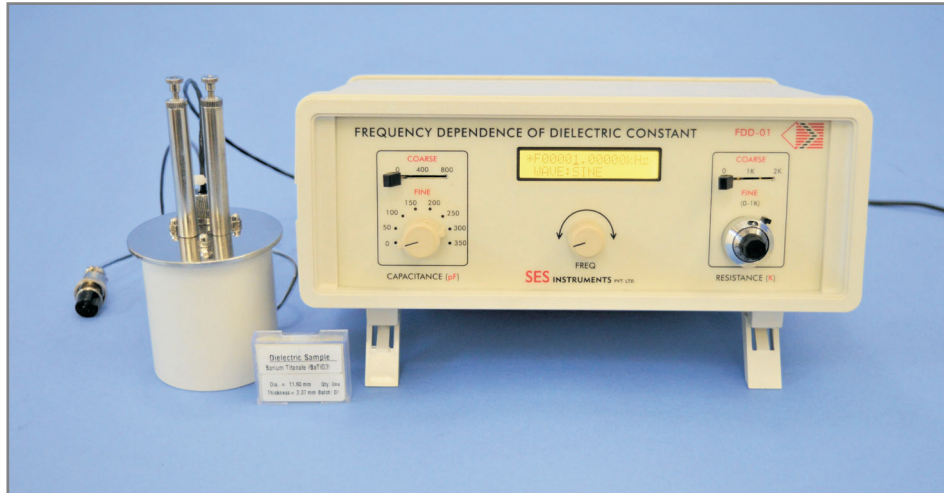

## Frequency Dependence of Dielectric Constant

- Capacitance measurement from 1-50kHz
- Dedicated Schering Bridge with built-in oscillator
- Computation of dielectric constant and loss factor
- Samples of Barium Titanate, MLCC and conventional capacitor
- Temperature variation studies with additional equipment oven and controller (optional)
- Essential accessory a CRO, for all the experiments



## Introduction

Barium Titanate,  $BaTiO_3$ , is a well known and widely investigated dielectric material. It is mainly used in capacitors due to its high dielectric constant. The dielectric properties of  $BaTiO_3$  are controlled by purity and microstructure which are dependent on the methods of preparation. The ability of the dielectric material to store energy is attributed to the polarization which can result in an increase in capacitance. It is well known that dielectric properties of every solid are very sensitive to the local electric field distribution in the sample. The temperature and frequency dependence of dielectric constant and loss can therefore give useful information about structure changes, transport mechanism and defect behaviour of a solid.

The present experiment is designed to study the variation of capacitance and dielectric loss as a function of frequency and temperature, to observe the change in dielectric constant

frequency dependent, the performance of a standard multi layer ceramic capacitor (MLCC) and conventional ceramic capacitors are not as prominent. A comparison of these three therefore is an interesting part of the study.

In addition, an optional experiment to study the variation of dielectric constant with temperature at different frequencies is also possible with the help of our PID controlled oven, PID-TZ

### Experiments

- Study of dielectric constant variation with frequency of a Barium Titanate sample
- Study of loss factor with frequency of a Barium Titanate sample
- Study of capacitance variation with frequency of an MLCC capacitor
- Study of capacitance variation with frequency of a conventional ceramic cap.

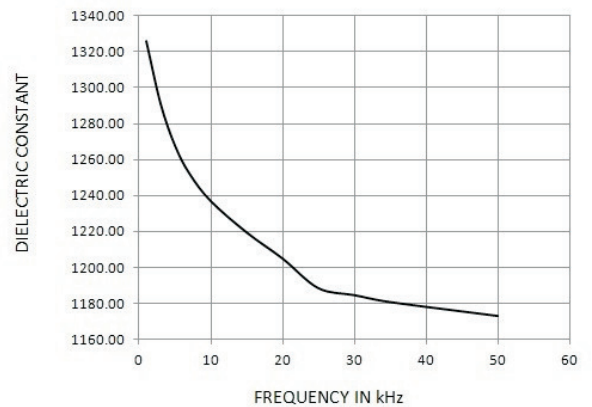


Fig 1  $\epsilon$  vs frequency of Barium Titanate sample

Typical result for Barium Titanate is shown in Fig 1.

### Optional Experiment (Using an additional PID-TZ)

- Study of Dielectric Constant of Barium Titanate as a function of temperature (curie point experiment) and its variation with frequency

Typical results are shown in Fig 2.

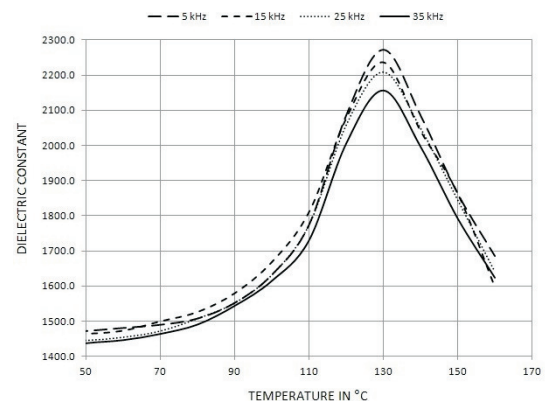


Fig 2  $\epsilon$  vs temperature at various frequencies of Barium Titanate sample

